
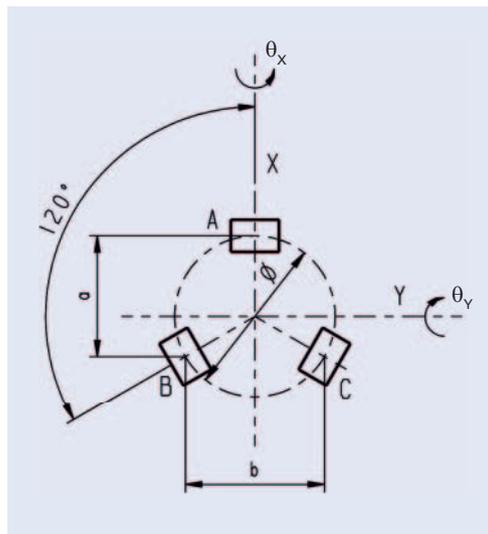


# Kinematics of Multi-Axis Tip/Tilt Systems

Piezo tip/tilt mirror systems from PI are based on parallel kinematics with a single movable platform for all directions of motion. The systems achieve a higher linearity than can be attained by switching two single-axis systems in succession, as is the case with galvanoscanners, for example, and therefore, are very compact.

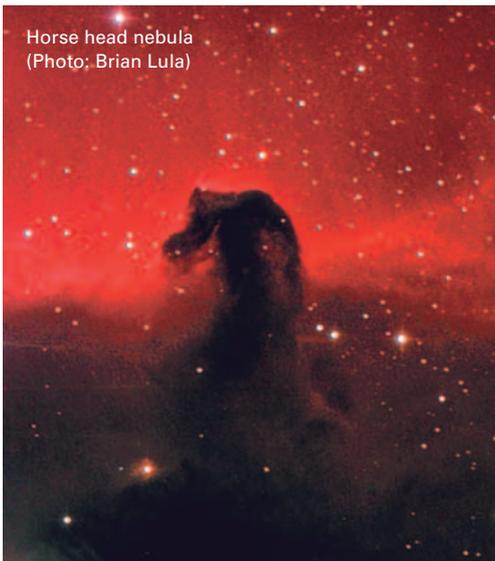

Arrangement of the actuators of a tripod piezo drive

The tilt angle and the travel in Z are calculated using the following formulas:

$$\theta_y = 2A - \frac{(B+C)}{2a}$$

$$\theta_x = \frac{(B-C)}{b}$$

$$Z = \frac{(A+B+C)}{3}$$

A, B, C is the linear displacement of the relevant actuators.

Piezo-actuated tip/tilt mirrors and platforms are suitable both for highly dynamic operation, such as tracking, scanning, image stabilization, elimination of drift and vibration, and for static positioning of optical systems and samples.

They allow for an optical beam deflection up to 100 mrad, extremely short response times from milliseconds to microseconds and resolutions down to nanoradians.

PI offers a large range from compact systems for laser beam steering up to large units used for astronomy.

### Tip/Tilt System with Tripod Piezo Drive

The platform is driven by three piezo actuators that are located in 120° angles to one another. By means of coordinate transformation, the motion can be split among the different actuators.


In addition to tilting, the platform may also be used linearly in Z direction, which is important, for example, for correcting optical path lengths (phase shifters).

### Tip/Tilt System with Differential Piezo Drive (Tetrapod)

The platform is driven by two pairs of piezo actuators located in 90° angles to one another. Four actuators are controlled differentially in pairs, depending on the tilt direction. The tilt axes  $\theta_x$  and  $\theta_y$  are arranged orthogonally so that a coordinate transformation is not necessary.

This concludes in an excellent stability in linear and angular positioning for a wide temperature range.

Just as the tripod, the differential version guarantees an optimum angular stability over a large temperature range. For position controlled versions, the differential evaluation of two sensors per axis provides an improved linearity and resolution.

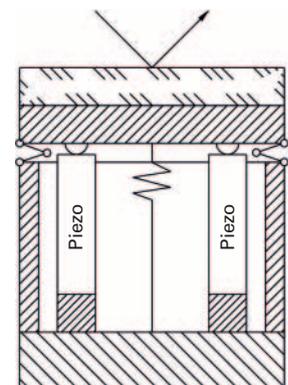

Principle of a tilt system with differential piezo drive

# Dynamics of a Piezo Tip/Tilt Mirror

The maximum operating frequency of a piezo tip/tilt system strongly depends on its mechanical resonant frequency. The properties of amplifier, controller and sensor are also important. To estimate the effective resonant frequency of the system – a combination of platform and mirror – it is necessary to calculate the moment of inertia of the mirror substrate first.

Moment of inertia of a rotationally symmetric mirror:

$$I_M = m \left[ \frac{2R^2 + H^2}{12} + \left( \frac{H}{2} + T \right)^2 \right]$$

Moment of inertia of a rectangular mirror:

$$I_M = m \left[ \frac{L^2 + H^2}{12} + \left( \frac{H}{2} + T \right)^2 \right]$$

with:

$m$  = mirror weight [g]

$I_M$  = moment of inertia of a mirror  
[g × mm<sup>2</sup>]

$L$  = mirror length orthogonally to tilt axis  
[mm]

$H$  = mirror thickness [mm]

$T$  = distance of pivot point to platform surface  
(see technical data of individual models)  
[mm]

$R$  = mirror radius [mm]

The resonant frequency of the system is calculated with resonant frequency of the platform (see technical data) and moment of inertia of the mirror substrate using the following formula:

$$f' = m \frac{f_0}{\sqrt{1 + I_M/I_0}}$$

Resonant frequency of a piezo tip/tilt system with mirror

with:

$f'$  = resonant frequency of platform with mirror [Hz]

$f_0$  = resonant frequency of platform without mirror [Hz]

$I_0$  = moment of inertia of platform (see technical data) [g × mm<sup>2</sup>]

$I_M$  = moment of inertia of mirror  
[g × mm<sup>2</sup>]



## Headquarters

### GERMANY

**Physik Instrumente (PI)  
GmbH & Co. KG**  
Auf der Roemerstrasse 1  
76228 Karlsruhe  
Phone +49 721 4846-0  
Fax +49 721 4846-1019  
info@pi.ws  
www.pi.ws

**PI miCos GmbH**  
Freiburger Strasse 30  
79427 Eschbach  
Phone +49 7634 5057-0  
Fax +49 7634 5057-99  
info@pimicos.com  
www.pi.ws

**PI Ceramic GmbH**  
Lindenstrasse  
07589 Lederhose  
Phone +49 36604 882-0  
Fax +49 36604 882-4109  
info@piceramic.com  
www.piceramic.com

© Physik Instrumente (PI) GmbH & Co. KG

All contents, including texts, graphics, data etc., as well as their layout, are subject to copyright and other protective laws. Any copying, modification or redistribution in whole or in parts is subject to a written permission of PI.

Although the information in this document has been compiled with the greatest care, errors cannot be ruled out completely. Therefore, we cannot guarantee for the information being complete, correct and up to date. Illustrations may differ from the original and are not binding. PI reserves the right to supplement or change the information provided without prior notice.

## Subsidiaries

### USA (East) & CANADA

**PI (Physik Instrumente) L.P.**  
Auburn, MA 01501  
www.pi-usa.us

### USA (San Francisco Bay Area)

**PI (Physik Instrumente) L.P.**  
Sausalito, CA 94965  
www.pi-usa.us

### ITALY

**Physik Instrumente (PI) S. r. l.**  
Bresso  
www.pionline.it

### FRANCE

**PI France SAS**  
Aix-en-Provence  
www.pi.ws

### JAPAN

**PI Japan Co., Ltd.**  
Tokyo  
www.pi-japan.jp

### CHINA

**Physik Instrumente  
(PI Shanghai) Co., Ltd.**  
Shanghai  
www.pi-china.cn

### SOUTHEAST ASIA

**PI (Physik Instrumente)  
Singapore LLP**  
Singapore  
www.pi-singapore.sg  
For ID / MY / PH / SG / TH / VNM

### KOREA

**PI Korea Ltd.**  
Seoul  
www.pikorea.co.kr

### USA (West) & MEXICO

**PI (Physik Instrumente) L.P.**  
Irvine, CA 92620  
www.pi-usa.us

### UK & IRELAND

**PI (Physik Instrumente) Ltd.**  
Cranfield, Bedford  
www.physikinstrumente.co.uk

### NETHERLANDS

**PI Benelux B.V.**  
Sint-Oedenrode  
www.pi.ws/benelux

### SPAIN

**Micos Iberia S.L.**  
Vilanova i la Geltrú  
www.pimicos.es

**PI Japan Co., Ltd.**  
Osaka  
www.pi-japan.jp

**Physik Instrumente  
(PI Shanghai) Co., Ltd.**  
Beijing  
www.pi-china.cn

### TAIWAN

**Physik Instrumente (PI)  
Taiwan Ltd.**  
Taipei  
www.pi-taiwan.com.tw